



BIOPHARMACEUTICAL DEVELOPMENT PROGRAM

SOP Title: Maintenance and Storage of Dissolved Oxygen and pH Probes
SOP Number: 12118
Revision: 05

TABLE OF CONTENTS

1. PURPOSE	1
2. SCOPE	1
3. RESPONSIBILITIES	1
4. MATERIALS AND REAGENTS.....	2
5. EVALUATION OF DISSOLVED OXYGEN SENSORS	2
6. PROCEDURE.....	3
7. STORAGE OF DISSOLVED OXYGEN SENSORS.....	4
8. EVALUATION OF PH PROBES	4
9. REPLACING A PH PROBE	4
10. STORAGE OF PH PROBES.....	4
11. DOCUMENTATION AND RECORDS.....	4
12. REFERENCES AND RELATED DOCUMENTS.....	5

1. PURPOSE

This Standard Operating Procedure (SOP) describes the maintenance and storage procedures for dissolved oxygen probes.

2. SCOPE

This SOP applies to Production personnel using reusable vessel mounted Dissolved Oxygen and pH Probes in BDP Upstream Production Areas.

3. RESPONSIBILITIES

3.1 Director, Late Process Sciences, Biopharmaceutical Development Program (BDP)

- Defines this procedure.

3.2 Manager, Large Scale Manufacturing, BDP

- Trains personnel in this procedure.

3.3 Area Supervisor, or designee(s)

- Implements this procedure.

BIOPHARMACEUTICAL DEVELOPMENT PROGRAM

SOP Title: Maintenance and Storage of Dissolved Oxygen and pH Probes

SOP Number: 12118

Revision: 05

3.4 BQA.

- Provides quality oversight of this procedure.

4. MATERIALS AND REAGENTS

Part Number	Description	BDP Approved Substitution Permitted?
21718	Mettler Toledo pO2 membrane kit (Small probes)	<input type="checkbox"/> YES <input checked="" type="checkbox"/> NO
20894	Broadley-James pO2 membrane kit	<input type="checkbox"/> YES <input checked="" type="checkbox"/> NO
21332	Broadley-James replacement cathode, pO2	<input type="checkbox"/> YES <input checked="" type="checkbox"/> NO
20554	Broadley-James pH probe replacement kit	<input type="checkbox"/> YES <input checked="" type="checkbox"/> NO
20632	pH Storage Solution	<input checked="" type="checkbox"/> YES <input type="checkbox"/> NO
NA	Oxy-probe polarizer (either cordless or corded style)	<input type="checkbox"/> YES <input checked="" type="checkbox"/> NO

5. EVALUATION OF DISSOLVED OXYGEN SENSORS

- 5.1 Prior to being placed into service for a run, perform a visual inspection.
- 5.2 Check for holes in the membrane tip and the condition of the outer cylinder O-ring.
- 5.3 The membrane must be replaced if there is a hole in it or media has migrated into the probe interior.
- 5.4 Failure of a probe to standardize or poor response and performance during a run also necessitates membrane replacement. It is recommended that for mammalian cell culture purposes, DO probes should have their membranes changed every 1-2 runs. The probe's response to testing prior to a run will help decide whether or not to replace the membrane.
- 5.5 The membrane must be replaced during product changeover.

SOP Title: Maintenance and Storage of Dissolved Oxygen and pH Probes
SOP Number: 12118
Revision: 05

6. PROCEDURE

6.1 Replacing Membranes

- 6.1.1 Unscrew the end cap from the sensor shaft and pull it from the electrode.
- 6.1.2 Pull the membrane cartridge off the anode/cathode assembly and discard membrane cartridge.
- 6.1.3 Rinse the anode/cathode assembly with WFI and dry it with a piece of tissue paper.

Note: Dark, irregular color of the silver anode inhibits probe function. If present, clean per manufacturer's instructions with a soft brush and rinse completely with WFI.

- 6.1.4 Replace all O-rings and elastomers.
- 6.1.5 Completely fill the membrane cartridge with polarographic oxygen electrolyte. Avoid touching the membrane. Gently tap the cap until air bubbles have been removed from the electrolyte.
- 6.1.6 With the membrane cartridge in a vertical position and the sensor positioned over the cap with the slot in the silver anode facing away from the operator, slip the membrane cartridge onto the anode/cathode assembly. Work any air away from the silicone sleeve. Allow air and excess electrolyte to escape from the anode groove. Continue pushing the membrane cartridge until it reaches the trapezoidal washer (if present).
- 6.1.7 Wipe away any expelled electrolyte solution from the electrode. Make sure that no electrolyte remains between the membrane cartridge and the trapezoidal washer.
- 6.1.8 Carefully slip the end cap over the fitted membrane cartridge and screw it into place.
- 6.1.9 Affix a label to the probe indicating "membrane changed", initials, and date.
- 6.1.10 Attach the polarizer and allow the probe to polarize for at least 2 hours before use.

BIOPHARMACEUTICAL DEVELOPMENT PROGRAM

SOP Title: Maintenance and Storage of Dissolved Oxygen and pH Probes
SOP Number: 12118
Revision: 05

7. STORAGE OF DISSOLVED OXYGEN SENSORS

- 7.1 When not in use, connect probes to a polarizer.
- 7.2 Store probes dry in such a fashion to protect the delicate membrane from damage.

8. EVALUATION OF PH PROBES

- 8.1 Prior to being placed into service for a run, perform a visual inspection.
- 8.2 Check for cracks in the glass that would compromise the ability of the probe to operate.
- 8.3 Failure of a probe to standardize or poor response and performance during a run also necessitates replacement of the probe.

9. REPLACING A PH PROBE

- 9.1 Unscrew the lock nut at the top of the probe housing.
- 9.2 Remove the pH probe from the housing and replace with an equal unit.
- 9.3 Replace all O-rings inside the housing, which are included in the pH probe replacement kit.

10. STORAGE OF PH PROBES

- 10.1 When not in use probes should be stored in a volume of pH storage solution to cover the exposed probe section.
- 10.2 pH storage solution should be changed on a monthly basis.

11. DOCUMENTATION AND RECORDS

- 11.1 Document cleaning on **Form 12149-01** as per **SOP 12149, General Cleaning of Process Equipment**.
- 11.2 Document maintenance such as membrane replacement in the fermentor/bioreactor logbook where the probe is used per **SOP 21531, Equipment/Facility Logs**.



BIOPHARMACEUTICAL DEVELOPMENT PROGRAM

SOP Title: Maintenance and Storage of Dissolved Oxygen and pH Probes

SOP Number: 12118

Revision: 05

12. REFERENCES AND RELATED DOCUMENTS

Document Number	Title
12149-01	Equipment Cleaning
12149	General Cleaning of Process Equipment
21531	Equipment/Facility Logs